

N605

4 Channel 200 MHz Constant Fraction Discriminator



Redesigning a classic: the new Constant Fraction Discriminator for fast high performance detectors

Features

- Four independent Constant Fraction Discriminators for fast detectors with sub-ns rise time
- Independently adjustable discriminators parameters (T, W, Z)
- Double pulse resolving time down to 4 ns
- Fine threshold adjustment from -15 to -1000 mV
- Fine walk correction with zero-crossing adjustment
- Exceptional low walk jitter < 50 ps
- Adjustable output signal width
- CFD monitor output
- Common veto input and individual gates for coincidence or anticoincidence options

Applications

- Fast Scintillators
- PMTs
- Silicon Detectors
- MCP

The N605 houses four independent 20% Constant Fraction Discriminators (CFD) in 1-unit wide NIM.

The module is intended for high resolution time measurements, where fast signals with rise time as low as 0.8 ns are involved.

Those may involve the use of radiation detectors made of fast scintillators, PMT, microchannel plates and fast semiconductor detectors.

The module guarantees exceptional low walk jitter, lower than 50 ps, for signals spreading over a wide range of amplitude, from -5 V to -50 mV.

The unit is provided with independent threshold discriminators to reject signal baseline noise. Front panel test point and screwdriver permits the precise measurement of the threshold, in the range from -15 mV down to -1 V. Each channel can be adjusted for the zero crossing level of the discriminator to minimize the time walk. Moreover the N605 has a CFD signal monitor output connector on the front panel to facilitate the correct adjustment of the working parameters.

Each channel has a fan-out of three timing output signals provided with NIM standard. The outputs can be selected to operate both in blocking and updating mode: the first is intended to minimize multiple triggering when slow scintillators are used, while the second is intended to reduce dead time in event of high rate condition. The board is provided with fast VETO input and individual GATE inputs to operate coincidence or anticoincidence logic conditions with the timing signal.



N605 rear view

Ordering Option

Code	Description
WN605XAAAAA	N605 - 4 Channel 200 MHz Constant Fraction Discriminator



News from Catalog web page
www.caen.it/news



n Small details
 Great differences



Copyright © CAEN SpA - 2015
 All rights reserved. Information in this publication supersedes all earlier versions. Specifications subject to change without notice.
 Printed in July 2015 - ADOCUME00116 - BF3268 - rev03

CAEN SpA
 Via Vetraia 11
 55049 - Viareggio • Italy
 Phone +39.0584.388.398
 Fax +39.0584.388.959
info@caen.it
www.caen.it

CAEN GmbH
 Klingenstrasse 108
 42651 - Solingen • Germany
 Phone +49.212.2544077
 Fax +49.212.2544079
info@caen-de.com
www.caen-de.com

CAEN Technologies, Inc.
 1140 Bay Street - Suite 2C
 Staten Island, NY 10305 • USA
 Phone +1.718.981.0401
 Fax +1.718.556.9185
info@caentechnologies.com
www.caentechnologies.com